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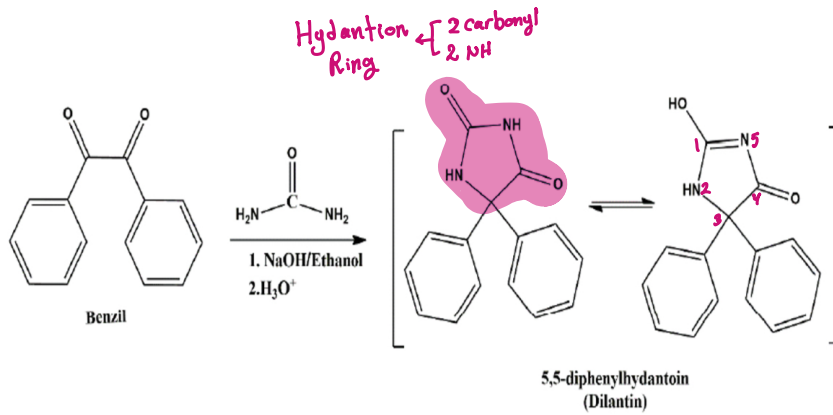
Step 3 : condensation of benzil with urea

- The condensation of benzil with urea, catalyzed by a strong base like sodium hydroxide (NaOH), is a reaction that synthesizes phenytoin (Dilantin), an anticonvulsant drug. The reaction involves the nucleophilic attack of deprotonated urea on one of benzil's carbonyl groups, followed by an intramolecular cyclization and a benzilic acid rearrangement that includes a-phenyl shift. The final product, phenytoin is typically isolated by acidifying the reaction mixture and collecting the resulting precipitate.

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
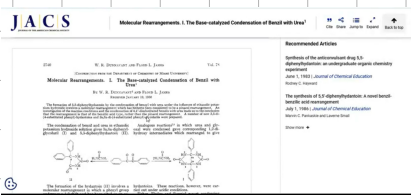
Step 3 : condensation of benzil with urea

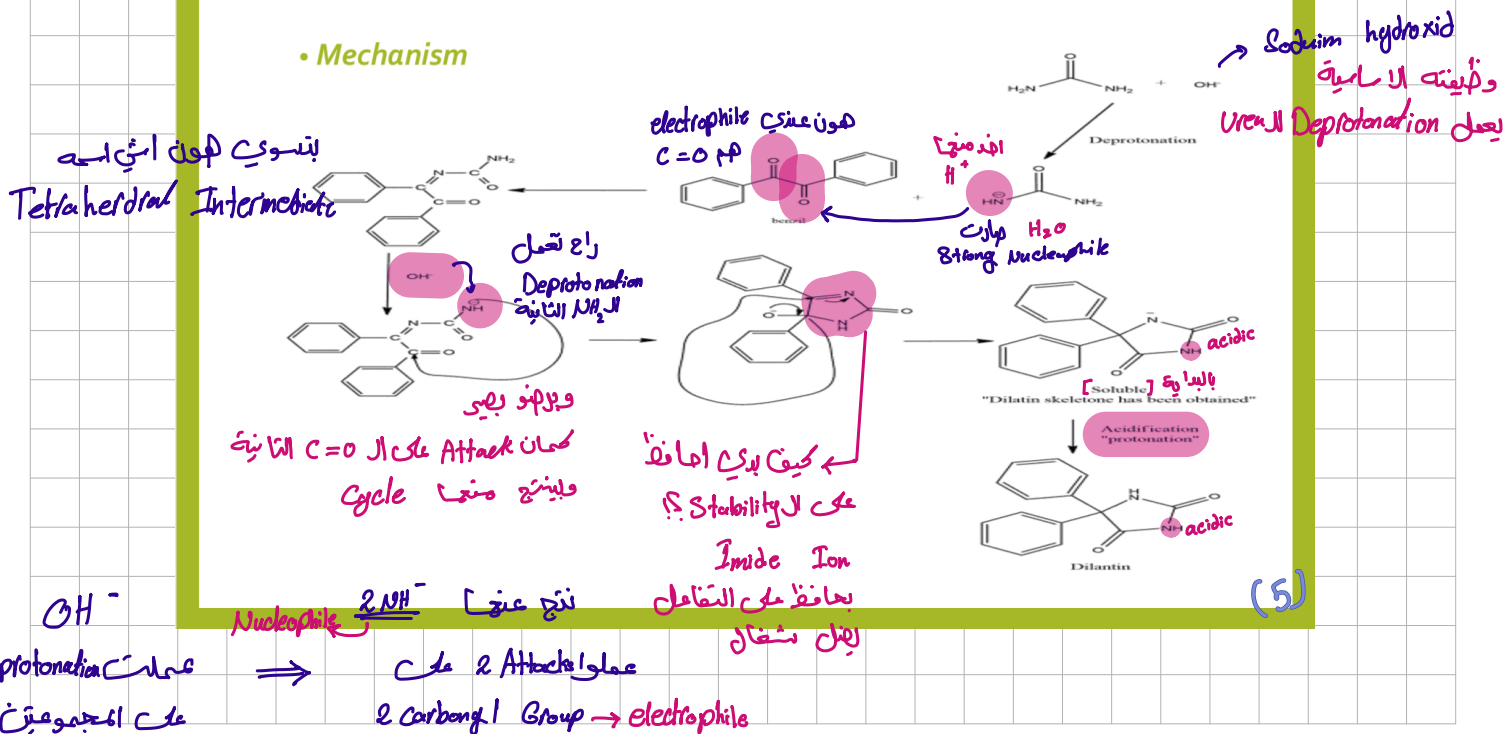• Mechanism

(5)

Phenytoin synthesis

Experiment 3 : Condensation of benzil and urea

① weigh 0.4 g of benzil compound prepared in the previous lab.

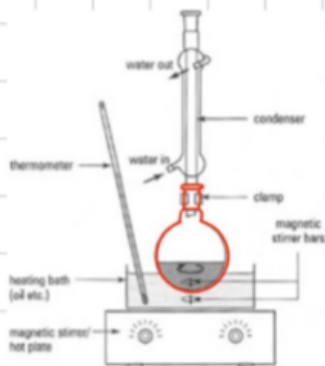
② In a round bottom flask add 0.4 g of the benzil compound, 200 mg urea, 6 ml absolute ethanol and shake, then add 1.2 ml of 30% aqueous sodium hydroxide and mix well.

③ attach the flask to the reflux condenser for 1 hour

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

Boiling
Chips

Reflux condenser

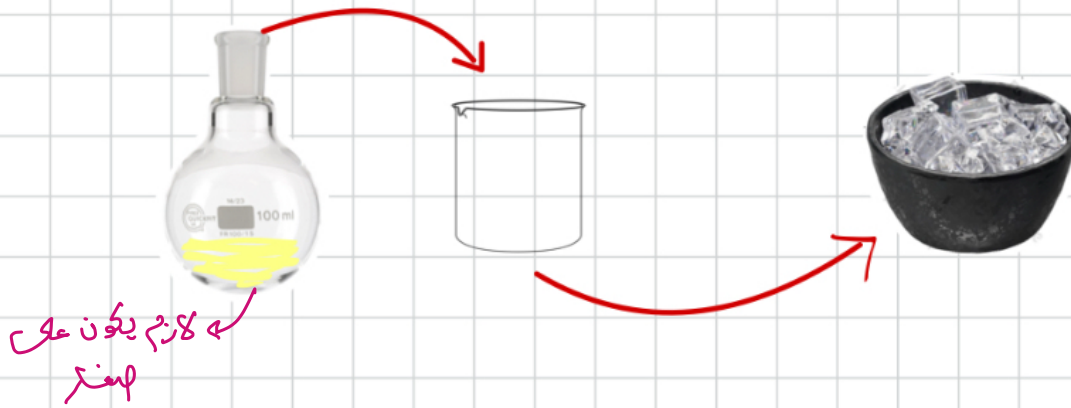③ Transfer your mixture to a 250 ml beaker and cool it in the ice bath .

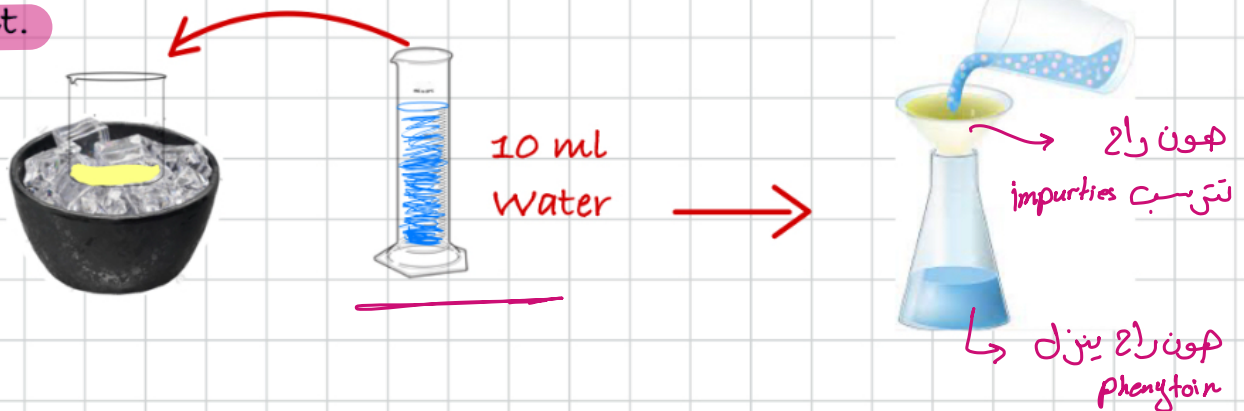④ Add 10 ml of water and filter the solution to remove the soluble side product.

⑤ Acidify your filtrate using 20% HCl to precipitate phenytoin (to get pH = 2 red on litmus paper)

⑥ cool it in the ice bath then collect your product using suction filtration (rinse with water)