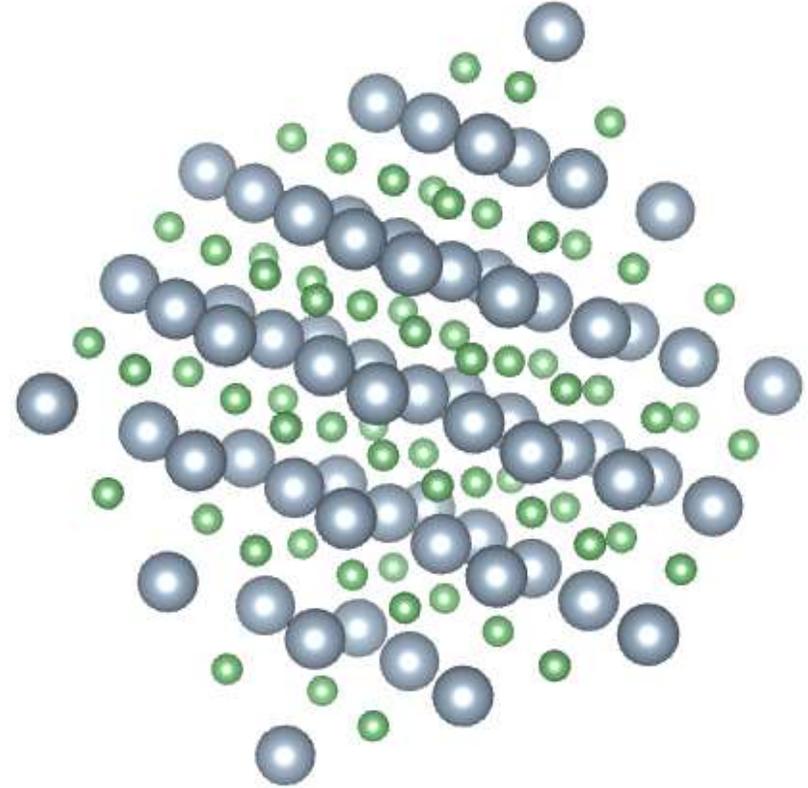
Melting Point Measurement (M.P.)

physical property

- ❖ Melting Point Definition.
- ❖ Reasons for Melting Point Measurement.
- ❖ Sample preparation.
- ❖ Consideration for M.P. measurement.

- Many organic compounds are solids at room temperature as a result of strong intermolecular forces which hold the individual molecules together in a crystal lattice.
- The nature and strength of these intermolecular forces are responsible for the observed differences in melting point.
- In general, if the forces are strong, the melting point will be high, and if they are relatively weak, the melting point will be low

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
Melting Point Definition.

- The temperature at which the first crystal just starts to melt until the temperature at which the last crystal just disappears. for example, M.P. 147-149°C. (from solid to liquid)

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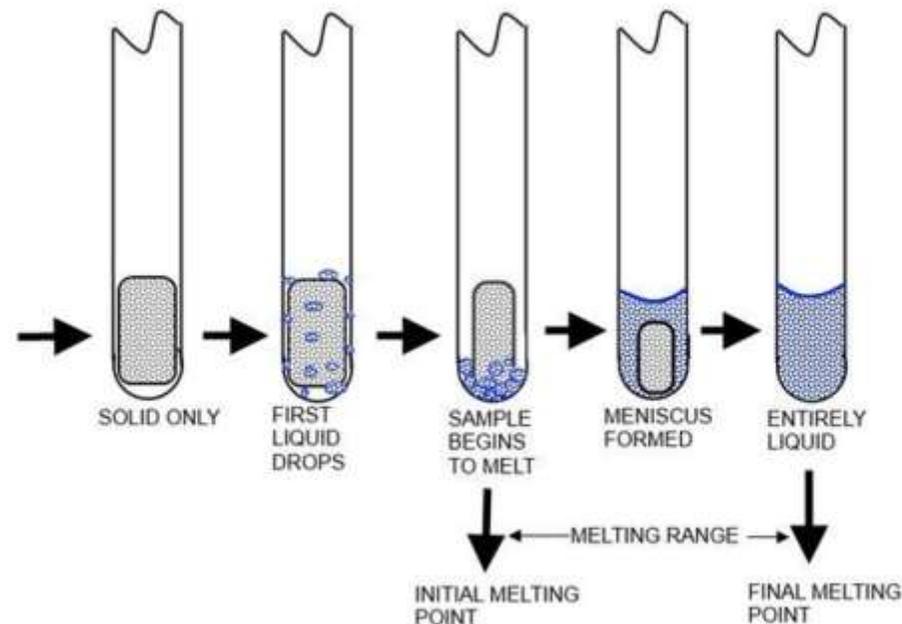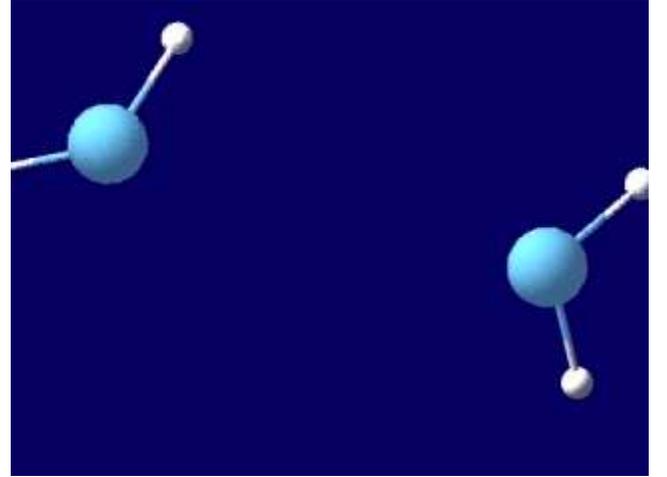
Soluble impurities affect the melting point of a solid in the following

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- a. **Lower the melting point of the substance**, with the upper limit considerably below the true melting point. The presence of an impurity in the molten compound, reduces its vapor pressure thus lowering the melting point of the compound. The greater the amount of impurity, the greater is the melting point depression.

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- b. **Broaden the melting point range**. Depending on the amount of impurity, the melting process may extend over a range of 2-20 oC or more.

- **Insoluble impurities** (*e.g., glass, sand ...etc.*) do not affect the melting point or the melting point range.

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Mixture melting points can be used in the following manner to determine whether two compounds are the same or different even though they have similar melting points.

EX

If a given organic compound (A) melts sharply at 120 °C,

and benzoic acid (compound B) also has a melting point of 120 °C. **Is compound (A) benzoic acid or a different compound?**

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- If **compound (A)** is benzoic acid, then a mixture melting point of (A) and (B) will melt sharply at 120 °C, i.e., the same as each individual compound alone.
- If, on the other hand, **compound (A) is not benzoic acid**, then the mixture melting point of (A) and (B) will be lowered and the melting range will be broadened. Since they are different compounds, each behaves as an impurity in the other.

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Reasons for Melting Point Measurement.

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1. Determination of **purity**. (pure substance has sharp range from 1 to 2

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a - measure then go to M.P. sheet

b - mix with Known substance with different ratio.

if two substances have the same M.P. then

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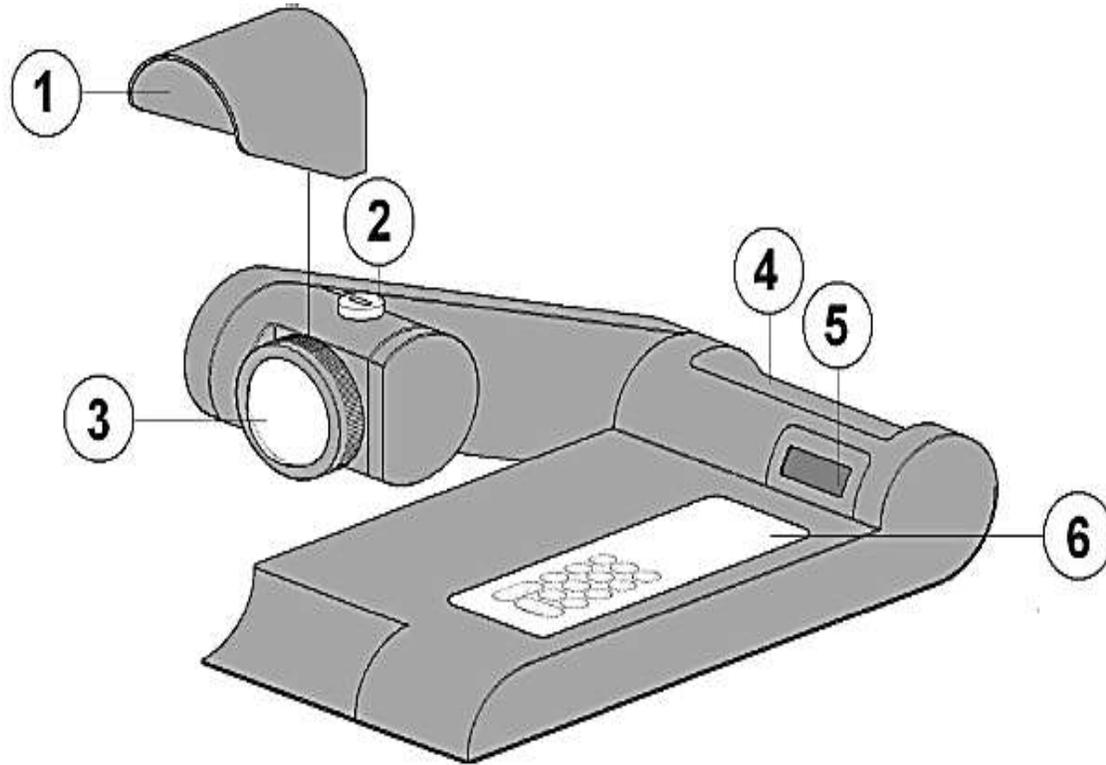
if not the same, you will notice lowering in the M.P. and Broad Range

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
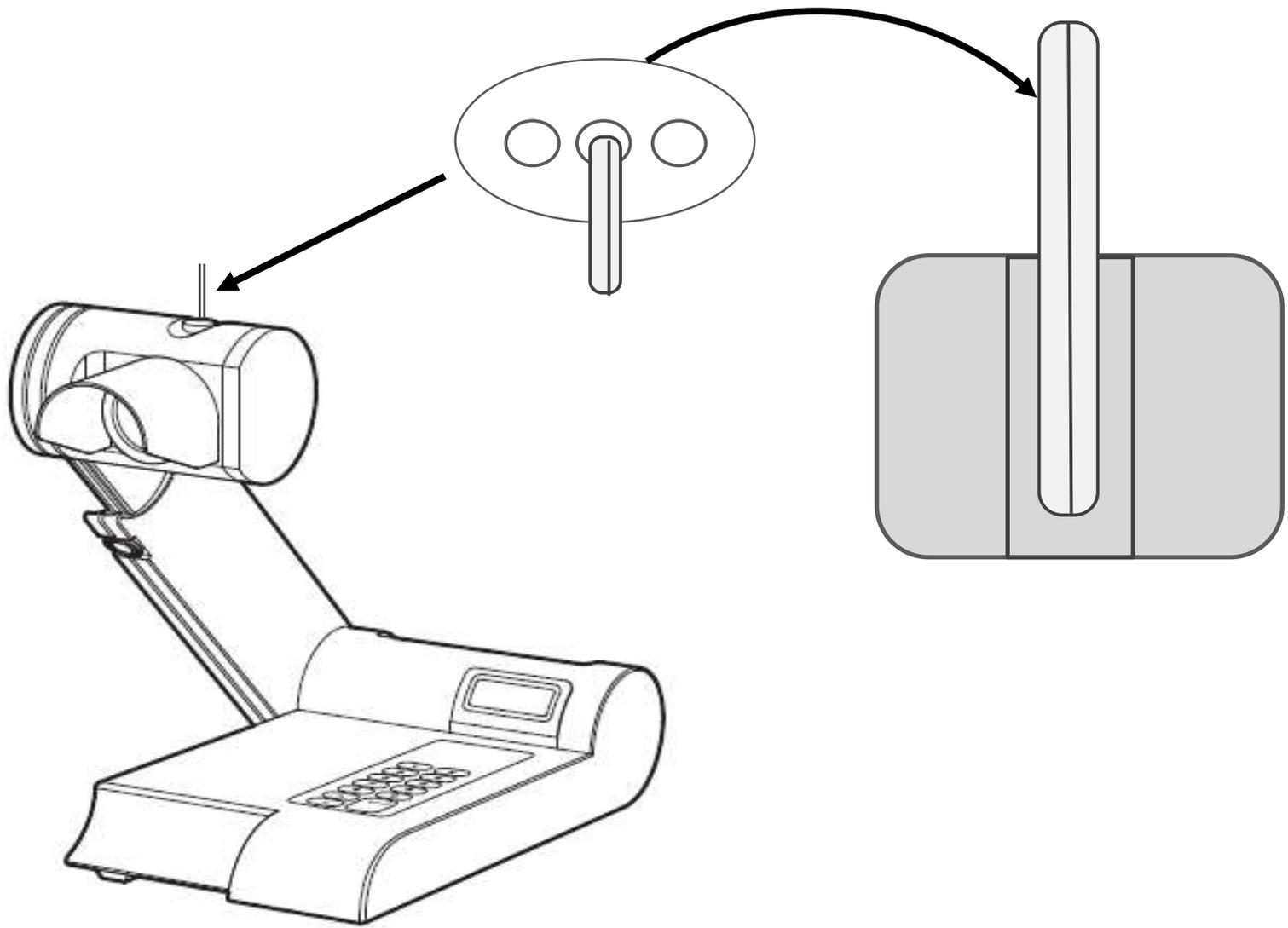
Sample preparation

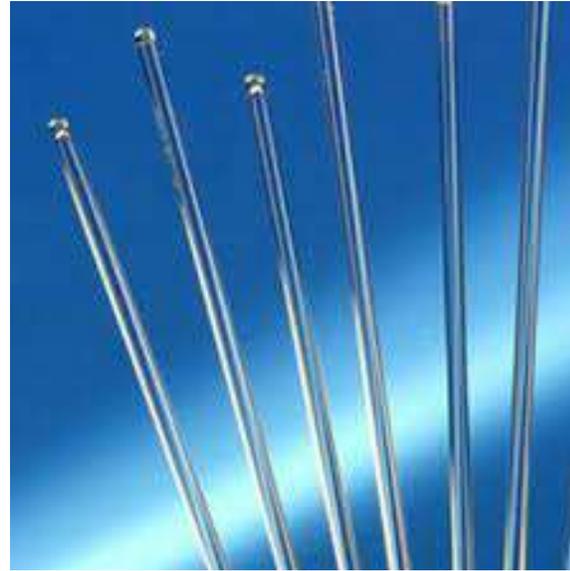GENERAL PROCEDURE

1. The digital melting point apparatus units must always be kept upright.
2. Place a small quantity (about 0.5 cm in tube) of the solid to be melted in a capillary tube (labeled melting point tubes).
3. Tap the closed end of the tube on the desk, clean the outside, and compact the solid down to the closed end of the melting point capillary tube.
4. Drop the tube (closed end down) down a section of glass tubing (see TA) to compact the solid in the bottom or closed end of the tube even more.
5. Place the tube loaded with the sample into the sample holder of the apparatus with the closed end down. The crystals can be ground up if they are too big to fit into the capillary tube.
6. Melting point capillary tubes are placed (closed end down) in the slots directly in front of the magnifying lens where they are viewed during melting. (Up to three samples can be viewed at once).

1. Record the temperature that the crystals begin to melt; crystals will look wet, (this is the melting start point), and the temperature at which the substance becomes a clear liquid; no solid material remaining (this is the melting end point).
2. Determine the melting range (Starting point – End point).
3. Calculate the melting point by taking the average point between the start point and the end point.

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The heating rate of the digital apparatus is adjusted by setting a temperature ramp along with a start and end temperature following the "Quick Start Instructions" on the front of the digital apparatus.

A ramp of 20 °C per minute will result in a rapid temperature rise while a ramp of say 2 °C per minute will give a slower rise that will more accurately measure the melting range of a solid.

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Procedure

Blueprint
EDUCATION
Quality • Precision • Focus

DETERMINATION OF MELTING POINT

Procedure

1. Powdering the compound and filling of capillary tube:
a) The thin walled capillary tube is sealed from one end by heating it on gas burner for a few seconds.

Blueprint Education © all rights reserved.

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Which of the melting point capillaries in the picture below is filled to the proper level for taking a melting point on a Mel-Temp melting point apparatus? (A, B)

A and B because we should put a small amount of the sample

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1. Use only the smallest amount that you can see melt. Larger samples will heat unevenly.

2. Pack down the material as much as you can. Left loose, the stuff will heat unevenly.

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3. Never remelt any sample. They may undergo nasty chemical changes such as oxidation, rearrangement and decomposition.

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4. Make up more than one sample. One is easy, two is easier. If something goes wrong with one, you have another. Duplicate, even triplicate runs are common.

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****Never repeat melting point's measurement with the same capillary as most of organic compounds decompose or react in other ways upon heating. This resulted in new materials with new melting points.**

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EXPERIMENTAL

MATERIALS NEEDED	<u>Equipment:</u> digital melting point apparatus <u>Glassware:</u> Capillary tubes (open one side only) <u>Chemicals:</u> pure and unpure mixures
-------------------------	--

- *DETERMINATION OF MELTING POINTS OF PURE COMPOUNDS*
- *IDENTIFICATION OF AN UNKNOWN*

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Measure the melting point by using the digital Melting point apparatus but you must first measure its melting range.

1. Tap a small amount of your unknown into two different capillary tubes. Just a **few crystals** are adequate. You may need to **grind some** of your unknown into a powder if it is too coarse to fit into the capillary tube.
2. Find the **melting point range of the pure unknown** substance by first quickly determining an approximate melting range **on a fast ramp (20 °C/min from 70-210 °C)**
3. Conduct a slow, careful **melting range with the second** capillary tube you prepared (use a ramp of 2 °C/min and start about 15 °C below the melting range to 10 °C above the range). Make sure the Digital Melting apparatus is below 70 °C before starting the first melting range and 10-20 °C below the compound's melting range before doing a slow careful melting range.

After determining the range, now you can calculate the approximate melting point midpoint

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DETERMINATION OF MELTING POINTS OF PURE COMPOUNDS

Each group should obtain a small amount (about 0.1 g) of one of the following solid compounds:

- | | | |
|------------------|-------------------|-----------------|
| 1. Cinnamic acid | 2. Salicylic acid | 3. Citric acid. |
| 4. Paracetamol | 5. Caffeine | 6. Urea |

Measure the melting point by using the digital Melting point apparatus but you must first measure its melting range.

1. Tap a small amount of your unknown into two different capillary tubes. Just a few crystals are adequate. You may need to grind some of your unknown into a powder if it is too coarse to fit into the capillary tube.
2. Find the melting point range of the pure unknown substance by first quickly determining an approximate melting range on a fast ramp (20 °C/min from 70-210 °C)
3. Conduct a slow, careful melting range with the second capillary tube you prepared (use a ramp of 2 °C/min and start about 15 °C below the melting range to 10 °C above the range). Make sure the Digital Melting apparatus is below 70 °C before starting the first melting range and 10-20 °C below the compound's melting range before doing a slow careful melting range.
4. After determining the range, now you can calculate the approximate melting point midpoint as follows:

$$\text{Melting Point Range} = (\text{Start Point} \rightarrow \text{End Point})$$

$$\text{Melting point Midpoint} = (\text{Start Point} + \text{End Point}) / 2$$

IDENTIFICATION OF AN UNKNOWN

Obtain an unknown (from instructor) and determine its melting point as described before. Using the melting points listed in the table on the previous page determine which possible compounds are within ± 10 °C of your unknown's melting range.

Table 1. Melting points of some organic compounds

Compound	mp (°C)	Compound	mp (°C)
Acetanilide	114	Maleic acid	135
Mandelic acid	117	Adipic acid	152
2-Naphthol	121	Citric acid	154
Benzoic acid	122	Salicylic acid	158
Urea	132	Benzanilide	161
Cinnamic acid	133	Sulfanilamide	165
Benzoin	133	p-Toluic acid	182